



## Gold Nanoparticles-NHS PRODUCT DATA SHEET

### Gold Nanoparticles-NHS

#### Description

Gold nanoparticles are widely used nanomaterials and generally referred to as colloidal gold in biological research. Colloidal gold markers generally has a particle size between 10 and 100 nm, and will show different colors with the change of particle size. Gold nanoparticles have excellent biocompatibility, rich surface modification properties, and unique optical properties related to the surfactant, shape, size, and structure of the nanoparticles. According to their different characteristics, it can be applied to various fields of biomedicine, such as medical testing, medical imaging, drug delivery, etc.

Abvigen provides a variety of gold nanoparticles, gold nanorods, gold nanocages, gold nanostars, gold nanobipyramids, and other products, the product particle size is optional, the concentration can be customized, the surface can be modified with different groups, and can be appropriately selected according to the customer's use.

For custom sizes, formulations or bulk quantities please contact our customer service department.

**Website:** [www.abvigen.com](http://www.abvigen.com) **Phone:** +1 929-202-3014 **Email:** [info@abvigenus.com](mailto:info@abvigenus.com)

#### Characteristics

Composition: Gold Nanoparticles-NHS

Shape: Spherical

Core diameter: 5 ~ 100 nm

Size: 1 mL

Amount: OD = 50

Surface: NHS

Buffer: DI Water

Form: Suspension



### Advantages

Results in covalently bound ligand and more stable conjugate.

Fast and convenient one-step conjugation reaction with no pre-activation requirements.

Spacer between the gold nanoparticle surface and conjugated ligand minimizes effects on the tertiary protein structure, which can lead to poor performing conjugates, a common problem seen in conjugates prepared by passive adsorption.

Precisely engineering surface coating that reduces non-specific protein binding in biological assays.

### Application

Ideal for development of protein gold conjugates for use in applications such as blotting, lateral flow assays, microscopy and transmission electron microscopy (TEM).

### Storage

This product should be stored at 4°C. **DO NOT FREEZE**. If stored unopened and as specified, Abvigen Gold Nanoparticles-NHS are stable for 3 months.

### Handling

When stored for a long period of time gold nanoparticles may sediment at the bottom of the vial, which is especially prominent for larger particle sizes. Prior to use, re-suspend the sedimented particles by thorough mixing until a homogenous solution is obtained.

### Note

These products are for R&D use only, not for drug, household, or other uses.

### NPS of Gold Nanoparticles, OD 50


Diameter	Peak SPR Wavelength	Optical density	Wt. conc	Size Dispersity %PDI	Particles/ml	Molarity mol/ml
5 nm	515-520 nm	OD 50	2.5 mg/ml	< 20%	1.98E+15	3.28E-09
10 nm	520 nm	OD 50	2.5 mg/ml	< 15%	2.47E+14	4.10E-10
20 nm	524 nm	OD 50	2.5 mg/ml	< 10%	3.09E+13	5.13E-11
30 nm	526 nm	OD 50	2.5 mg/ml	< 6%	9.15E+12	1.52E-11
40 nm	530 nm	OD 50	2.5 mg/ml	< 4%	3.86E+12	6.41E-12
50 nm	535 nm	OD 50	2.5 mg/ml	< 4%	1.98E+12	3.28E-12

60 nm	540 nm	OD 50	2.5 mg/ml	< 4%	1.14E+12	1.90E-12
70 nm	548 nm	OD 50	2.5 mg/ml	< 4%	7.21E+11	1.20E-12
80 nm	553 nm	OD 50	2.5 mg/ml	< 4%	4.83E+11	8.02E-13
90 nm	564 nm	OD 50	2.5 mg/ml	< 4%	3.39E+11	5.63E-13
100 nm	572 nm	OD 50	2.5 mg/ml	< 4%	2.47E+11	4.10E-13

#### Gold Nanoparticles Centrifugation Parameters

Particle Size	Speed (g)	Time (min)
5 nm	100000	30
10 nm	17000	60 (~ 50% recovery)
20 nm	6500	30
30 nm	4500	30
40 nm	2500	30
50 nm	2000	30
60 nm	1125	30
80 nm	400	30
100 nm	400	30

#### TEM of Abvigen gold nanoparticles of different size

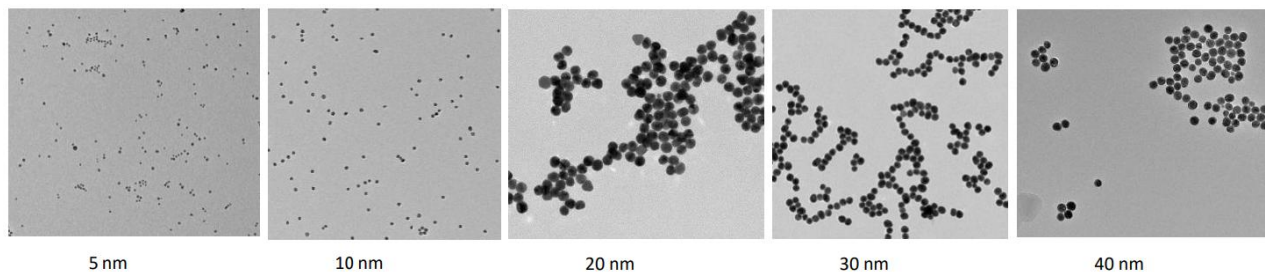

#### Ordering Information

Website: [www.abvigen.com](http://www.abvigen.com)

Phone: +1 929-202-3014

Email: [info@abvigenus.com](mailto:info@abvigenus.com)